OPSIS LiquidLINE KjelROC Analyzer

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THE OPSIS LIQUIDLINE LABCONNECT SOFTWARE

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ADAPTIVE WORKFLOW AND GLP REQUIREMENTS

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OPsis LIQUIDLINE MAINTENANCE PROGRAM

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KJELROC ANALYZER, KD-310-C1/C2

Analytical Data	
Titration System	Colorimetric Titration. Unique Predictive Titration system automatically adjust titration speed on the actual distance from the end point. Steam stirrer function ensure safe distillation to avoid dangerous exothermic reactions. PTFE tubing used throughout to eliminate formation of air bubbles between burette-titration vessel.
Sampler System	It is possible to install the KjelROC Analyzer with an optional Sampler system
Measuring Range w/o burette refill	0,1 - 225 mg N
Recovery	> 99.5% (1-200 mg N)
Burette volume	Interchangeable burette 50 ml (automatic refill during analysis)
Dosing accuracy	1.95 µl/step
Accuracy/Precision %rsd	0.75% at N levels 1 - 225 mg
Distillation Speed	40 ml/min at 230V
Start time for Analysis	Instant start of analysis, no delay between analyses
Distillation Time	3.5 - 4 min (10 mgN), 5-5.5 min (225 mgN)
Technical Data	
Operating Temperature	5°C - 40°C, max 80 % relative humidity
Power Supply	190-240 VAC, 50-60 Hz, 10A
Power consumption	max 2200 W
Cooling Water consumption	1,75 l/min at 20° C
Steamgen. water consumption	approx. 200 ml/analysis
Dimensions (WxHxD) and Weight	430 x 700 x 330 mm, 30 kg
Graphics Display	7" 800x480 color touchscreen
Safety	Safety sensors (Tube in place-, Protection- and Service door), Steam generator sensors, Cooling water Measurement, Drip tray, Tank level detectors, Maintenance warning systems
Electronic Steam Regulation	10-100%
Distillation Head	Easily exchangeable Splashhead with only one tube connector made of glass for high precision (B1) or Polypropylene for durability (B2).
Password Protection	Three Levels with separate Managers menu
Included Software	OPSIS Labconnect Standard software (can later be upgraded to LabConnect LIMS). OPSIS LiquidLINE Transfer Utility supplied for iPhone and iPad. Well documented open transfer protocols and example code for Excel spreadsheet provided.
Communication	Ethernet (RJ-45) interface. Included KjelROC router for wireless, 802.11 b/g.
Reagent Level Warning Data	4 x RJ-11 (external tank levels), 1 x RJ-11 (titration tank), OPSIS LiquidLINE flexible level detectors (optional accessory)
Automatic Distillation and Titration	Programs for automatic control of cooling water, dilution water, sodium hydroxide and receiver solution as well as automatic emptying of tube and titration vessel. Automatic control of water cooling, selectable delay times and cleaning program.
Programs	Up to 64 customized programs
Manual Control	Possibility to bypass automatic titration system to allow for manual titration.
Daily Maintenance	Manual and automatic cleaning programs

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Specifications subject to change without notice

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